

ASEPTIQUIK® S SERIES CONNECTORS

AseptiQuik® S Connectors provide quick and easy sterile connections for small-flow applications, even in non-sterile environments. The "FLIP-CLICK-PULL" design of AseptiQuik S enables users to easily transfer small volumes of media with less risk of operator error than with traditional methods. The connector's genderless and robust design provides reliable performance without the need for clamps, fixtures or tube welders. Biopharmaceutical manufacturers can now make 1/8", 1/4" and 3/8" hose barb and 1/4" and 3/4" sanitary sterile connections with the quality and market availability they expect from the leader in single-use connection technology.

SPECIFICATIONS

OPERATING PRESSURE

Up to 60 psi, 4.1 bar

OPERATING TEMPERATURE

39°F to 104°F (4°C to 40°C)

STERILIZATION

- ☐ **Gamma:** Up to 50kGy irradiation
- ☐ **AutoClave High Temp (HT) Version:** Up to 266°F (130°C) for 60 minutes

TERMINATIONS

1/8", 1/4" and 3/8" ID hose barb (3.2mm, 6.4mm and 9.5mm), 1/4" and 3/4" sanitary and MPC insert

MATERIALS

Main Components:

Polycarbonate (white)

Pull Tabs/Caps:

Polycarbonate (blue, standard version)

Polycarbonate (white, HT version)

Seals:

Silicone (clear), platinum-cured

Membrane:

Polyethylene (standard version)

Hydrophobic polyethersulfone (HT version),

PTFE strip sticker

WARNING: Pressure, temperature, chemicals, and operating environment can affect the performance of couplings. It is the customer's responsibility to test the suitability of CPC's products in their own application conditions.

FEATURES


- Genderless →
- FLIP-CLICK-PULL →
- CPC Click →
- AQS-MPC Combination →
- AQS ¼" Sani with Smooth Bore →

BENEFITS

- Eases single-use systems specifications with one-part number for both halves
- Intuitive three-step connection process reduces risk of operator error
- Audible confirmation of connection with no additional hardware required
- Enables the ability to change a BPC or other single-use system with open format connections to closed systems
- Minimizes transitional flow disruptions throughout upstream processing

TYPICAL FLOW RATE:

Cv Value Range: 0.19 - 1.74
for AseptiQuik S

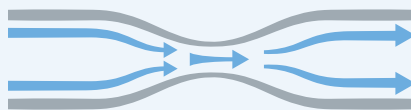

Cv values represent the approximate expected flow rate in gallons per minute of water at room temperature for a 1 PSI pressure drop. The flow is generally constrained by the smallest diameter, which in some cases will be the termination diameter and not the Nominal Flow Path.

NOTE

Validation and Extractables data can be requested at cpcworldwide.com/AseptiQuik-S

DID YOU KNOW

Did you know that the AseptiQuik S is perfect for simplifying the process to pull flask samples from your bioreactor?

Scan code to visit webpage



SCAN

cpcworldwide.com/AseptiQuik-S



COLDER PRODUCTS COMPANY
U.S.A.
PHONE: +1 (651) 645-0091
FAX: +1 (651) 645-5404
TOLL FREE: +1 (800) 519-7633

COLDER PRODUCTS COMPANY GMBH
Germany
PHONE: +49 6026 9973-0

COLDER PRODUCTS COMPANY LIMITED
ShangHai, China
PHONE: +86 21 2411 2666
TOLL FREE: +86 400 990 1978

ASEPTIQUIK S SERIES DIMENSIONS

POLYCARBONATE with blue pull tabs - *For gamma irradiation applications.*




TERMINATION	METRIC EQ.	PART NO.	A	B	C
1/8" HOSE BARB	3.2 mm	AQS17002	2.25" (57.2 mm)	1.30" (33.0 mm)	0.50" (12.7 mm)
1/4" HOSE BARB	6.4 mm	AQS17004	2.25" (57.2 mm)	1.45" (36.8 mm)	0.65" (16.5 mm)
1/4" SANITARY	6.4 mm	AQS33004	2.25" (57.2 mm)	1.50" (38.1 mm)	0.70" (17.8 mm)
3/8" HOSE BARB	9.5 mm	AQS17006	2.25" (57.2 mm)	1.45" (36.8 mm)	0.65" (16.5 mm)
3/4" SANITARY	19.1 mm	AQS33012	2.25" (57.2 mm)	1.60" (40.6 mm)	0.80" (20.3 mm)
MPC INSERT		AQS17MPC	2.25" (57.2 mm)	1.49" (37.9 mm)	0.69" (17.5 mm)

POLYCARBONATE HT with white pull tabs - *For autoclave or gamma irradiation applications.*



TERMINATION	METRIC EQ.	PART NO.	A	B	C
1/8" HOSE BARB	3.2 mm	AQS17002HT	2.25" (57.2 mm)	1.30" (33.0 mm)	0.50" (12.7 mm)
1/4" HOSE BARB	6.4 mm	AQS17004HT	2.25" (57.2 mm)	1.45" (36.8 mm)	0.65" (16.5 mm)
1/4" SANITARY	6.4 mm	AQS33004HT	2.25" (57.2 mm)	1.50" (38.1 mm)	0.70" (17.8 mm)
3/8" HOSE BARB	9.5 mm	AQS17006HT	2.25" (57.2 mm)	1.45" (36.8 mm)	0.65" (16.5 mm)
3/4" SANITARY	19.1 mm	AQS33012HT	2.25" (57.2 mm)	1.60" (40.6 mm)	0.80" (20.3 mm)
MPC INSERT		AQS17MPCHT	2.25" (57.2 mm)	1.49" (37.9 mm)	0.69" (17.5 mm)

PRODUCT DIMENSIONS

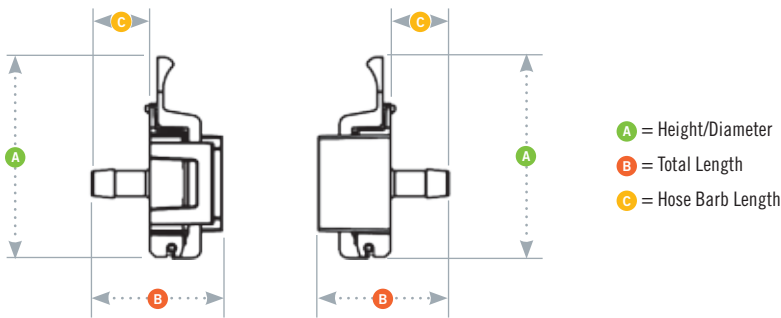

NOTES

WARRANTY: All sales are subject to Colder Products Company's limited express warranty set forth in the CPC catalog. Contact your local distributor or CPC Customer Service for warranty provisions.

Warning: Due to the wide variety of possible fluid media and operating conditions, unintended consequences may result from the use of this product, all of which are beyond the control of CPC. It is the user's responsibility to carefully determine and test for compatibility for use with their application. All such risks shall be assumed by the buyer.

COPYRIGHT © 2021 BY COLDER PRODUCTS COMPANY. COLDER PRODUCTS COMPANY, COLDER PRODUCTS AND CPC ARE REGISTERED

TRADEMARKS WITH THE US PATENT & TRADEMARK OFFICE